LOVIBOND®
MINIKIT

GB: Organo-Phosphonate
D: Organo-Phosphit
F: Organo-Phosphite
I: Organo-fofrito

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GB

Introduction

The Organo-Phosphonate Minikit provides a simple means of monitoring organophosphonate levels over the range 1 - 20 mg/l and is based on a titration. The test has been developed for use with commercially available organo-phosphonate products based on amino trimethylphosphonic acid and hydroxyethane diphosphonic acid- the two compounds most commonly used in these formulations.

Contents

titrator tube

direct reading titrator

Organo- Phosphonate No.1 tablets, 50 pieces in foil

Organo- Phosphonate No.2 solution, 100 ml

operating manual in 8 languages

Instructions

Range 1 - 20 mg/l

Sample Size 10 ml

1. Take a sample of the water under test in the titrator tube, filling to the 10 ml mark.
2. Add one Organo-Phosphonate No. 1 tablet, and shake the tube gently to disintegrate the tablet.
3. Fill the direct reading titrator with Organo-Phosphonate No.2 solution (Fig. 1).
Ensure there are no air bubbles trapped in the titrator. Adjust so that the solution level coincides with the zero mark (Fig.2).
4. Insert the titrator through the hole in the cap of the titrator tube (Fig.3).
5. Gently depress the plunger on the titrator, adding the solution a drop at a time and shaking the tube gently to ensure mixing.
6. Continue adding the solution until the colour changes
from green to purple.
7. Observe the mark corresponding to the solution in the titrator (Fig. 4). This figure represents the active organo-phosphonate concentration in the sample as milligrams per litre.

Cleaning

Thoroughly rinse out the titrator tube and the direct reading titrator after each use.

Important

1. The results of this test are expressed in terms of mg/l (ppm) active organo-phosphonate. Commercially available products are normally sold as aqueous formulations with a given active organo-phosphonate content. To calculate the dosage of commercial products required from the test results, regard must be made of the active content of the product in use.
2. For most accurate results it is recommended that a blank be carried out on the water without organo-phosphonate compound, or on the makeup water to the system and the result of this test subtracted from the test reading obtained with the cooling water.
Substances reported to interfere with this test are aluminium, polyphosphate, chelating agents and iron in high concentrations.

Use of the direct reading titrator

Fig. 1 Fill the titrator with reagent solution

Fig. 2 Adjust reagent level to zero mark

Fig. 3 Insert titrator into tube and add reagent, one drop at a time

Fig. 4 Read result from titrator mark

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solution d'Organo-Phosphonate No.2, 100 ml

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